

Coffee, Sugar and NanoCuvette™ One

NanoCuvette™ One offers a quick and easy way of detecting sugar in your coffee.



1. Coffee

Coffee is a darkly coloured drink, brewed on roasted coffee beans from the Coffea plant. It originates from tropical Africa, but is now cultivated in several equatorial countries across South-America, Asia and Africa, and is one of the most popular drinks in the world. Coffee can be prepared and served in a variety of ways. It is most often served warm, but may also be served iced. Some people take it with milk or cream, others with sugar.

2. Sugar detection

Sugar content is typically determined based on refractive index measurements. For pure solutions of a known compound the refractive index measurement can be used to determine the exact concentration of the specific sugar. For mixed solutions with high sugar content, such as coffee with sugar, the refractive index measurement can be used to detect the sugar and give an approximate measure of the sugar content

3. Principle

Conventionally a specialized refractometer has been required for the refractometric measure of sugar. However, with the innovative NanoCuvette™ One, a nanosensor is installed in a cuvette, allowing measures of sugar to be carried out in a standard spectrophotometer. Because sugar does not absorb light in the UV-visible spectrum, traditionally, spectrophotometers have been unable to measure sugar in coffee.

4. Safety precautions

This method does not entail any safety precautions. Please refer to common laboratory practices.

5. Measurement

5A. Materials and apparatus

The only apparatus required is a standard spectrophotometer and a computer with internet access. For each measurement a NanoCuvette™ One is required.

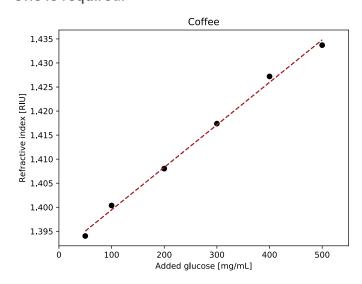


Figure 1 The sensor response, shown as refractive index, is linearly proportional to sugar concentration, which the free software will display.

5B. Sample preparation

Coffee is prepared according to standard methods, and 1 to 500 mg/mL sugar is added.

www.nanocuvette.com Page 1

Application Note

5C. Measurement procedure

The spectrophotometer and computer are switched on and the free NanoCuvette™ One software is opened. For each sample a new NanoCuvette™ One is used according to the software guidance. Following the last measurement the software will present the data as a variable of absorbance or refractive index, depending on the experimenter's preferred choice of representation (Figure 1).

6. Contact

Copenhagen Nanosystems Diplomvej 381 DK-2800 Kgs. Lyngby

Tel: +45 36 99 27 46 info@cphnano.com www.nanocuvette.com

Updated September 2017



Figure 2 NanoCuvetteTM One measures refractive index and absorbance on the same sample, using only a standard spectrophotometer and the free NanoCuvetteTM One software.

www.nanocuvette.com Page 2